WATER BATH VERIFICATION PROCEDURE #22

December 1, 1996

Equipment Checked: WATER BATHS (AASHTO T 209 and T 245) (ASTM D 3666)

Purpose:

This method provides instructions for checking water temperatures in a bath.

<u>Inspection Equipment Required:</u>

1. Thermometer

Tolerance:

```
25 degrees C ± 1 degree C
60 degrees C ± 1 degree C
```

Procedure:

Place thermometer into bath and read temperature. Adjust as necessary.

EQUIPMENT CALIBRATION RECORD

Calibration Procedure No. 22 Water Bath (AASHTO T 209)

Date:	Calibrated by:
Previous Calibration Date:	Next Due
Frequency: 6 Months	
Calibration Equipment	Serial Number
Thermometer	
Heater Setting	
Thermometer Reading	
Comments:	

5/9/2002